Miltenyi Biotec

gentleMACS™ Octo Dissociator with Heaters

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ANCELL

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Figure 4.1: Front of the gentleMACS Octo Dissociator with Heaters.

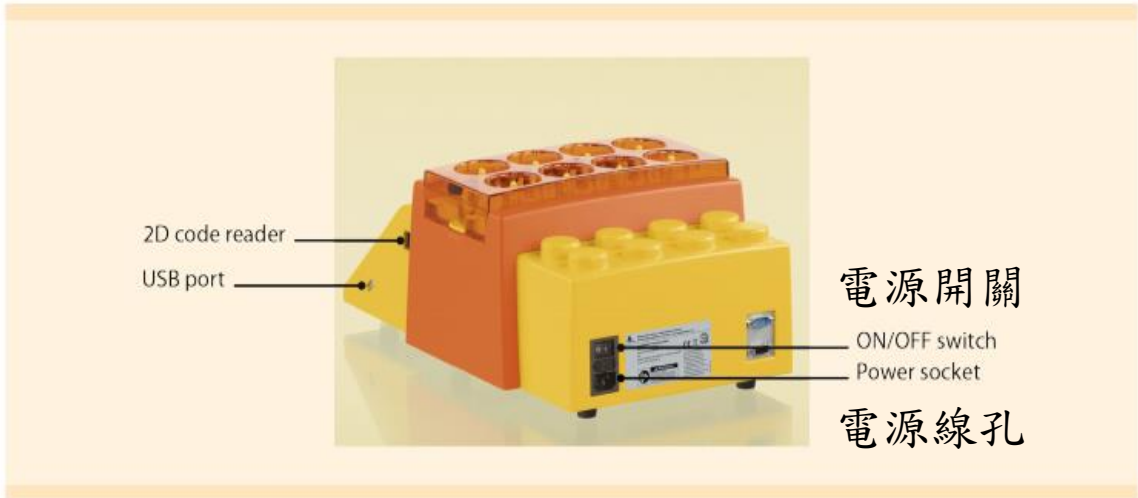
Figure 4.2: Back of the gentleMACS Octo Dissociator with Heaters.

Figure 4.3: Back of the gentleMACS Octo Dissociator with Heaters, Safety Shield removed.

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Icon	Folder	Description
	Miltenyi	The Miltenyi folder contains gentleMACS Programs provided by Miltenyi Biotec. The lock on the Miltenyi folder indicates that programs in this folder cannot be edited or deleted.
	Templates	The Templates folder contains gentleMACS Programs provided by Miltenyi Biotec. Programs in this folder can be edited but must be renamed and saved in a User folder.
	User - 1	Five User folders are available to manage user-defined programs.
	User - 2	
	User - 3	
	User - 4	
	User - 5	
	Favorites	Links of programs in the Favorites folder enable quick access.
	USB stick	A USB stick can be installed to transfer programs to and from a folder.

Table 4.2: Overview of folders and respective icons.

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Figure 4.4: Main menu of the gentleMACS Octo Dissociator with Heaters.

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